

Lynx EVO - SmartCam5/50


IMPORTANT: Before installing software, make sure you read the instructions and follow Lynx EVO assembly steps in the User's Guide.

- 1 Unpack
- 2 Assemble Lynx EVO
- 3 Install Software


Min. Recommended System Specs:

Windows 10 (32/64-bit), Intel® Core i5 - 2.5 GHz, 4GB RAM / 200 GB HD

1. Uninstall any previous version of IDS uEye, Pylon and ViCapture software from your PC.
2. Do not plug SmartCam 50 camera into the PC
3. **Install IDS Peak** – “ids_peak_2.1.0.0.exe”
 - Follow On-Screen Wizard
 - Select Setup type as Custom Install

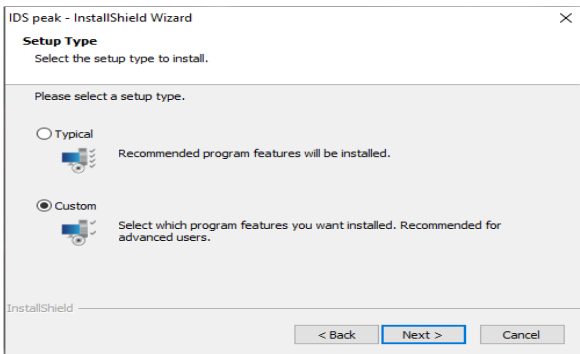

- At “Select Features” window, select the option “**Support for the uEye cameras**”

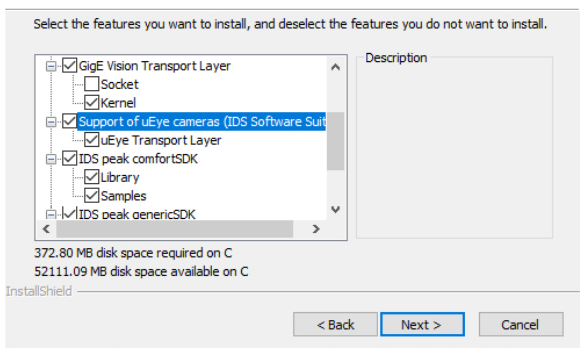

- Complete the installation process and “**Restart the PC**”

4. **Install uEye** v4.951 or later – “uEye_49510.exe”
 - Complete the installation process and “**Restart the PC**”
5. Connect the SmartCam 5 /50 camera to USB 3.0 or SS port on PC using the provided USB cable
6. Lynx EVO with SmartCam 5/50 is now ready for use.
7. Open **IDS Peak Cockpit** from your desktop